TiO₂ Nanotube Synthesis

Fluoride etching of Ti metal substrate to form TiO2 nanotubes

The procedure described in detail below is taken from the following publications. If this procedure is followed, please cite appropriately.

- Meekins, B. H.; Kamat, P. V., Got TiO2 Nanotubes? Lithium Ion Intercalation can Boost Their Photoelectrochemical Performance Three-Fold. *ACS Nano* **2009**, 3, 3437–3446. NDRL 4821 http://dx.doi.org/10.1021/nn900897r
- Baker, D. R.; Kamat, P. V., Disassembly, Reassembly and Photoelectrochemistry of Etched TiO2 Nanotubes. *J. Phys. Chem. C* **2009**, 113, 17967-17972. NDRL 4816 http://dx.doi.org/10.1021/jp9065357
- Baker, D. R.; Kamat, P. V., Photosensitization of TiO2 Nanostructures with CdS Quantum Dots. Particulate versus Tubular Support Architectures. *Adv. Funct. Mater.* **2009**, 19, 805-811. NDRL 4771 http://dx.doi.org/10.1002/adfm.200801173
- 1. Sonicate Ti foil (0.8cm x 4.0cm) in 2-propanol for 1 hour, store in acetone until use
- 2. Fill a 3-arm electrochemical cell with electrolyte solution
 - Electrolyte Solution (200mL batch)
 - i. Dissolve 0.6686g NH₄F in 4mL H₂O
 - ii. Add solution to 196mL ethylene glycol
- 3. Place Ti metal foil (cathode) and a platinum counter electrode (anode) in electrolyte solution.
- 4. Slowly increase bias voltage between electrodes from 0V to 60V (about 1 V/s ramp)
 - Bubbles should begin to form off of counter electrode
- 5. Allow etching to take place for desired amount of time (generally 1-2 hours depending on desired nanotube length)
- 6. Turn off power supply and remove etched Ti foil. Rinse with EtOH
- 7. Sonicate in EtOH for 2 seconds (further sonication leads to detachment of the nanotubes from the Ti substrate
- 8. Dry in air and sinter at 450°C for 3hrs